

# Installation Instructions for Micro Series Mixer Cartridges (binary & ternary housings)

1. To replace an existing mixer cartridge, disassemble the mixer housing using 5/8" & 1" wrenches (or use two adjustable wrenches).
2. Store the old cartridge in a manner that will protect the end seals from being scratched or damaged. Also, make sure that no particles are allowed to block the tiny holes in the ends of the cartridge.
3. Insert the replacement cartridge and reassemble the housing. Please note that there may be a white residue on the housing threads. This is an inert Teflon coating that prevents galling or seizing of the threads, and should not be removed.
4. Tighten the housing finger tight until housing makes contact with cartridge, then tighten an additional **1/4 turn** with wrench. **DO NOT TIGHTEN FURTHER OR MIXER PERFORMANCE MAY BE REDUCED.**

